

**0.9 % NaCl in tap water**

9 gm NaCl /L tap water

or

216 gm NaCl / 24 L tap water

**Volumes:**

**Total: 24 L** in the labeled carboy in Burrill 230A

\*\*\*\*\*

**2.0 % Urea in tap water**

20 gm Urea / L tap water

or

480 gm Urea / 24 L tap water

**Volumes:**

**Total: 12 L** in the labeled carboy in Burrill 230A

\*\*\*\*\*

For each Section you will need:

4 Fleakers each with 1 L 2% urea and 1.7 grams flavoring

4 Fleakers each with 1 L 0.9% NaCl and 1.7 grams flavoring

4 Fleakers each with 1 L Tap water and 1.7 grams flavoring

4 Fleakers each with 1 L Tap Water

Put in Cooler in 217 BH for at least 3 hours before lab